WiseStir[®] & ***WiseTherm***[®]
Digital Hotplate / Hotplate Stirrer
/ Magnetic Stirrer

Written for
MSH-20D, MSH-30D, MS-20D, HP-20D & HP-30D

Operation Manual
(Version 1.3.2)

DAIHAN Scientific Co., Ltd.

1

Introduction

Thank you for purchase of "**Digital**" *WiseTherm*[®] Hotplate, *WiseStir*[®] Magnetic Stirrer & Hotplate Stirrer.

Those products have superior features as follows;

- Mixing liquids, Stirring, Heating and Controlling
- Digital Feedback control by High-Performance Microprocessor
- Superior Temperature Uniformity: Less than 3% on the plate at current temperature (*HP-20D, HP-30D, MSH-20D & MSH-30D*)
- The Ergonomic Article for precise-control of HEAT, STIRRING -speed, and TIME Jog-Shuttle switch (Turn + Push)
- Accurate Temp. Control of the sample in the vessel using Optional PT100 Sensor (*HP-20D, HP-30D, MSH-20D & MSH-30D*)
- User's self-compensation function: to control Difference between Real Temp. and Indicated Temp. ($\pm 10^{\circ}\text{C}$) (*HP-20D, HP-30D, MSH-20D & MSH-30D*)
- Ceramic Coated Plate: Chemical/Acid Resistance
- Smart Design of Body: made by Aluminum Die-Casting (*MS-20D, MSH-20D & HP-20D*)
- Storage function for the set values: Temperature, RPM and Timer
- Digital LCD with Back-Light Function
- Digital Timer function included: 99hr 59min settable
- Alarm function: for Error Status of Timer-end
- Locking Mode supported for Experimental Safety (Input to Jog-Shuttle can be disabled)
- Free Voltage (AC 80V~240V, 50Hz ~ 60Hz) by SMPS (*MS-20D*)

Package

Contents

2

MSH-20D Digital Hotplate Magnetic Stirrer	<input type="checkbox"/>
MSH-30D Digital Hotplate Magnetic Stirrer	<input type="checkbox"/>
MS-20D Digital Magnetic Stirrer	<input type="checkbox"/>
HP-20D Digital Hotplate	<input type="checkbox"/>
HP-30D Digital Hotplate	<input type="checkbox"/>
SS100, Direct Contact Temperature Sensor (PT100)	<input type="checkbox"/>
SS110, Direct Contact Temperature Sensor (PT100)	<input type="checkbox"/>
SS200, Direct Contact Temperature Sensor, Teflon Coated (PT100)	<input type="checkbox"/>
SS210, Direct Contact Temperature Sensor, Teflon Coated (PT100)	<input type="checkbox"/>
RD100, Stand Rod	<input type="checkbox"/>
RD200, Stand Rod	<input type="checkbox"/>
CL 220, Holder + Clamp for the Temp. Sensor	<input type="checkbox"/>
Power Code	<input type="checkbox"/>
Operation Manual	<input type="checkbox"/>

3

Appearance

Each unit of *WiseTherm*[®] Hotplate, *WiseStir*[®] Magnetic Stirrer & Hotplate Stirrer has an independent model. The appearance of *WiseStir*[®] Hotplate Magnetic Stirrer, *MSH-20D* is shown in <Figure 1>.

<Figure 1> Appearance of *MSH-20D*

Models

- (a) *HP-20D*, *HP-30D* : Digital Hotplate
- (b) *MS-20D* : Digital Magnetic Stirrer
- (c) *MSH-20D*, *MSH-30D* : Digital Hotplate Stirrer

Operations

4


Multifunctional Digital Feedback Controller

<Figure 2> Controller of *WiseStir*[®] Digital Hotplate Stirrer

- ① Operation Status on LCD Display
- ② Indication of Set Value
- ③ Indication of Current Value
- ④ Power Bar Graph: indicating the power of motor operating
- ⑤ LED indicating Heating Operation. **(MS-20D n/a)**
- ⑥ LED indicating Stirring Operation. **(HP-20D, HP-30D n/a)**
- ⑦ LED indicating Timer Operation.
- ⑧ Mode Switch
- ⑨ Jog Shuttle Switch: Turn + Push

Back plane of *HP-20D*, *HP-30D*, *MS-20D*, *MSH-20D* & *MSH-30D*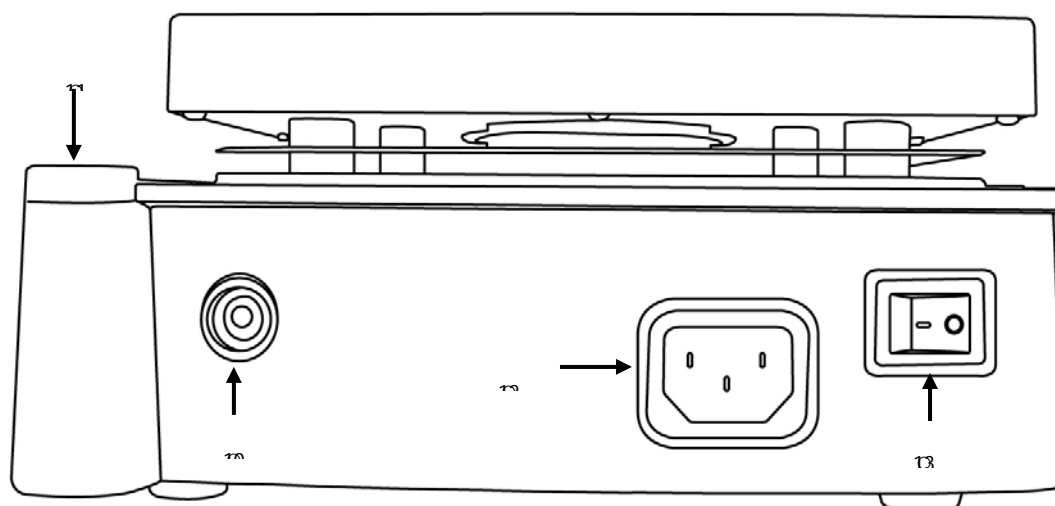

10 Socket for connecting SS100, SS110, SS200 or SS210 (Optional External PT100

Sensor) (**MS-20D n/a**)


11 Holder for the Support Rod, RD100 or RD200

12 Power Code connector

13 Power Switch

Operation guide

1. Power-on: Turn on the Power Switch (ⓓ), and then get into its default mode.
2. Mode Selection: You can change the mode by click MODE switch (⑧).

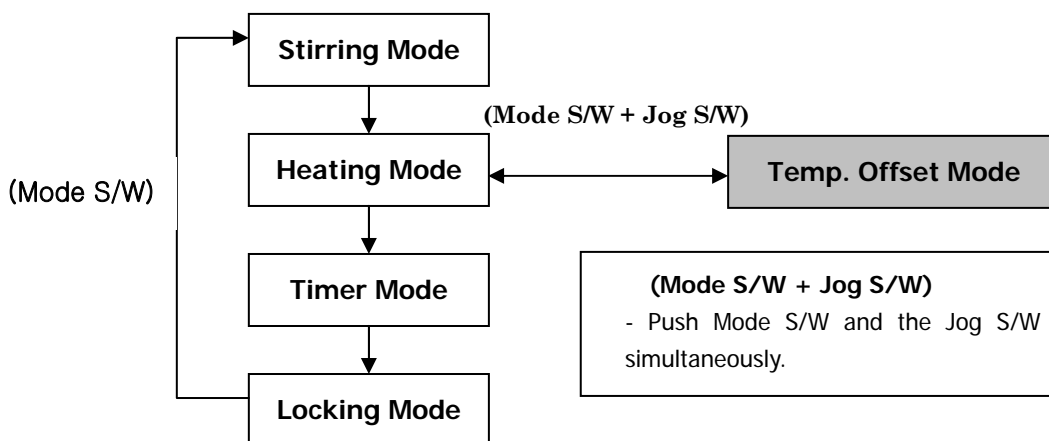We support two different temperature control system as follows whether usage of the external sensor, SS200 (optional PT100 sensor) or SS2100 (optional Teflon coated PT100 sensor) or not. **(MS-20D n/a)**

- A. If you don't connect PT100 sensor, the temperature is controlled by the measured value of hotplate itself.
- B. If you connect PT100 sensor, the temperature is controlled by digital feedback and the current value of temperature (③) shows the actual value of solution.

The external PT100 sensor is recognized automatically when you plug it in the device.

Set Value and Current Value indicate different value at each operation mode, and it is summarized on the Table 1.

Setting Mode	Set Value	Current Value	Control target
Stirring Mode	Set(Saved) Speed (rpm)	Current Speed (rpm)	Stirring Speed (rpm)
Heating Mode	Set(Saved) Temperature	Current Temperature	Temperature of Hotplate or Sample
Off Set	Compensated Temperature	Off set	Calibration of temperature

Timer Mode	Set(Saved) Time	Remained Time	Operation time
Locking Mode	Current Temperature	Current Speed	Display only

<Table 1> The meaning of digit at the each mode

3. Stirring Mode / Heating Mode

- A. After turning on the power, select one between Stirring Mode and Heating Mode and you can confirm the LED of selected mode is flickering. On the panel, the Set Value shows zero or non-zero and the Current Value shows zero if the device is not operated.
 - i. If the set value is non-zero, it means someone has saved the shown value previously. You can use the value simply by clicking (pushing) the jog shuttle at the each mode. Or, you can change the value by turning the jog shuttle at the each mode and click the knob. Then, the device begins to revolve at the set speed or to heat to be reached at the set temperature according to the mode.
 - ii. If the set value is zero, it means any values were not saved and you can adjust the set RPM or temperature by turning the jog shuttle. Then, push the knob and the device begins to revolve at the set speed or to heat to be reached at the set temperature according to the mode.
- B. RPM (Temperature) adjustment: When you turn the jog shuttle for adjustment of RPM or temperature, the set value is change firstly and the current value goes after the set value. Then, the current value is reached to the set value, and it means that *WiseTherm*[®]/*WiseStir*[®] Hotplate Magnetic Stirrer series are feedback controlled by digital controller accurately. (Control resolution: 5 rpm for speed, 0.5°C for temperature)
- C. Pause and Restart Function: If you click the jog shuttle during stirring operation in RPM mode, stirring is temporary stopped (Pause Status). If you click again the knob, the device is restarted the operation.
- D. Save the current set value: You could save the current RPM (temperature) in your operation by clicking the jog shuttle for one second. This value is saved in the device while it is turned off and you can reuse it without setting values.

4. Timer Mode

- A. You can use the internal digital timer to preserve your current operation status during the set time. After termination of the set time, the operation is automatically stopped. So you have to set up the operation modes (the requested RPM and temperature) before you set up the timer.
- B. Set the Timer: To set up the timer, you have to enter the Timer Mode by clicking the mode button, and then turning the jog shuttle to adjust timer.

Click (push and then release immediately) the jog shuttle after turning to set required value. When you enter the Timer Mode, you could see a time set before. You are able to use that value itself, or to modify by turning the jog-shuttle.

- C. Start and pause timer: Click the jog-shuttle to start the timer. During the operation of timer, the LCD shows the remained time. Click again to pause the timer.
- D. Restart overall operations: After termination of set time, the operation is automatically paused. And, you need to release the pause of each operation mode by pushing the jog shuttle if you want to restart overall timer operations.

5. Locking Mode

After completion of set up the RPM, temperature and timer, you can set the Locking (Safe) Mode by clicking the mode button. In the Locking Mode, the LCD shows the current temperature and RPM in measuring value digit. In this mode, all inputs are disabled using the jog shuttle. If you want to release this mode, click the mode select button.

6. Temperature Offset (MS-20D n/a)

Offset Function is addition or subtraction of indicated temperature. It is very convenient function to synchronize displayed temperature of the plate and sample with the measured value by the User.

- A. Press Mode Switch and Jog Shuttle simultaneously in the Heating Mode (Mode S/W + Jog S/W).
- B. In this status, it is changed into "Offset" Mode and the user can adjust the value by pushing the knob once, then the offset would be setup.
- C. After the setup, the temperature on LCD display indicates the compensated temperature.
- D. At this time, press the mode switch and jog shuttle simultaneously, then escape from Offset Mode.

7. Storage of the set value

WiseTherm[®]/*WiseStir*[®] Digital Hotplate and Magnetic Stirrer have a "Memory" function to save each set value. The user can save the set values to be used frequently and does not have to set the same value every moment.

- A. Set the value of each mode (except Locking mode) using Jog Shuttle.

Press the knob over one second, then the set value is blinking on LCD three times and it means that the value has been saved.

8. Power off

After you complete to use the device, you must turn off the power switch for your safety (⚠).

Safety

Instructions

5

Do not modify any part of the machine - the mechanism part or electrical part. Modification made by non-technical person may cause a serious problem of the machine.

Do not use flammable material.
Flammable materials or Chemicals must not be used in order to avoid an explosion or a fire.

The white plate could be raised up to 350°C during heating, and touching this surface may cause a serious injury.
If you need to touch it, you should have sufficient cooling time down to

70°C on the LCD display before touching.

The recommended power supply is AC 100-230V, 50/60Hz, Single phase.

The machine must be standing on a stable, even surface that is as sleep-free as possible.

Avoid allowing objects to push or strike the hotplate or stirring plate of the machine.

You have to use Hotplate magnetic stirrer at indoor environment, to preserve temperature between 5°C and 40°C.

The hotplate is made of ceramic-coated aluminum. Although the ceramic coating prevent the plate from erode by acid chemical, pay attention not to be stained with acid or basic material.

Please be careful that liquid should not penetrate into the part of control panel or back plane.

Please use the machine in clean environment for use with long lifetime.

You must not wrap the plate in aluminum foil, which can be damage the machine critically because of overheating of the body.

※ All circuits and electronic parts used for this product are DAIHAN's own technology, so please do not make it repaired by anyone who isn't our own staff or technician authorized by DAIHAN.

6

Specifications

WiseTherm[®] HP-D Digital Hotplate

		HP-20D	HP-30D
Plate Size (mm)		180 X 180	260 X 260
Temp.	Range (°C)	Max. 380	
	Accuracy	±0.3°C at Set Temperature	
Temp. Control Resolution (°C)		0.5	
Heating Power (W)		600	1200
Heating Power (W)		Specially Designed Heating Module (High Uniformity: Temp. Difference 3% less)	
Controller		Digital Feedback Controller with Jog Shuttle Switch (Turn + Push)	
Display		Digital LCD with Back-Light Function	
Timer		99hr 59min (Continuous function included)	
Material	Plate	Ceramic Coated (Chemical / Acid Resistance)	
	Body	Powder Coated Aluminum	Powder Coated Steel
Dimension (w X d X h mm)		206 X 307 X 99	286 X 387 X 99
Weight (Kg)		2.8	3.2
Others		Storage Function (Temp., Timer), Locking Mode (Jog-Shuttle Input disabled) User Self-compensation function Alarm (Error Status and Timer-end)	
Power Supply*		AC 110 V, 60 Hz or AC 230 V, 50 / 60 Hz	

WiseStir[®] MS-D Digital Magnetic Stirrer

		MS-20D
Stirring Capacity (H₂O)		Max 20Lit.
Motor Rating (W)		9 / 4
Speed Range (rpm)		80 ~ 1,500
Speed Control Resolution (rpm)		5
Speed Accuracy (%)		± 5
Controller		Digital Feedback Controller with Jog Shuttle Switch (Turn + Push)
Display		Digital LCD with Back-Light Function
Timer		99hr 59min (Continuous function included)
Plate Size (mm)		180 X 180
Material	Plate	Ceramic Coated Aluminum (Chemical / Acid Resistance)
	Body	Powder Coated Aluminum
Dimension (w X d X h mm)		206 X 307 X 99
Weight (Kg)		3.1
Others		Storage Function (RPM, Timer) Locking Mode (Jog-Shuttle Input disabled) Alarm (Error Status and Timer-end)
Power Supply*		Free Voltage (AC 80 ~ 240 V, 50 / 60 Hz)

WiseStir[®] MSH-D Digital Hotplate Stirrer

		MSH-20D	MSH-30D
Plate Size (mm)		180 X 180	260 X 260
Stirring Capacity (H₂O)		Max 20Lit.	
Motor Rating (W)		9 / 4	
Speed Range (rpm)		80 ~ 1,500	
Speed Control Resolution (rpm)		5	
Temp.	Range (°C)	Max. 380	
	Accuracy	±0.3°C at Set Temperature	
Temp. Display Resolution (°C)		0.1	
Temp. Control Resolution (°C)		0.5	
Heating Power (W)		600	1200
Heating Plate		Specially Designed Heating Module (High Uniformity : Temp. Difference 3% less)	
Controller		Digital Feedback Controller with Jog Shuttle Switch (Turn + Push)	
Display		Digital LCD Display with Back-Light Function	
Timer		99hr 59min (Continuous function included)	
Material	Plate	Ceramic Coated (Chemical / Acid Resistance)	
	Body	Powder Coated Aluminum	Powder Coated Steel
Dimension (w X d X h mm)		206 X 307 X 99	286 X 387 X 99
Weight (Kg)		3.3	3.5
Others		Storage Function (Temp., RPM, Timer) Locking Mode (Jog-Shuttle Input disabled) User Self-compensation function Alarm (Error Status and Timer-end)	
Power Supply*		AC 110 V, 60 Hz or AC 230 V, 50 / 60 Hz	

CERTIFICATE

of

Quality Tested by Manufacturer and Free Warranty up to 2 Years

For DAIHAN Laboratory Equipment

Limited Warranty

1. This instrument has gone through our testing process about all functions before shipment.
2. This instrument is warranted for two years from purchasing date.
3. A breakdown caused by customer's misuse or natural disaster is excepted from this warranty.
4. This certificate is only for international customers.

Item	
Model No.	
Serial No.	
Date	
Customer	

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